


A microscopic view of cells, likely lymphocytes, with a large cell in the upper right and several smaller cells scattered across the field. The background is a soft, out-of-focus mix of pink, purple, and blue.

Life Sciences
Cytokine assays



cisbio

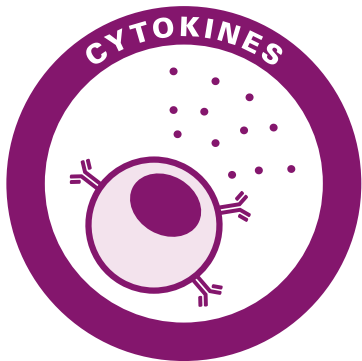

A full range of cytokines

Cisbio's kits cover the full range of cytokines: interleukins, interferons (IFN), chemokines, TNF and lymphokines.

All our assays are cell based and measure cytokine release in cell supernatant.

They have been validated in several cell lines, primary cells and PBMCs to ensure relevance, and their low sample requirement enables the quantification of multiple cytokines in parallel from the same supernatant.

The HTRF® protocol makes them a fast alternative to ELISA or other measurement methods for researchers in the fields of oncology, inflammatory diseases or any other inflammation process.

This leaflet presents the analytical performances of each assay, the associated standard curve* and the list of available references.**

* To learn more about data processing and 4PL $1/y^2$ fitting, we invite you to visit our dedicated webpage at <http://www.cisbio.com/4PL-regression>


** Bulk size kits are available - please contact us

STRONG ASSAYS


- ✓ Lyophilized standards
- ✓ Low sample volume: 16 μL
- ✓ Same buffers
- ✓ Single protocol
- ✓ 4PL $1/y^2$ fitting

RELY ON VALIDATION


Supernatant from:

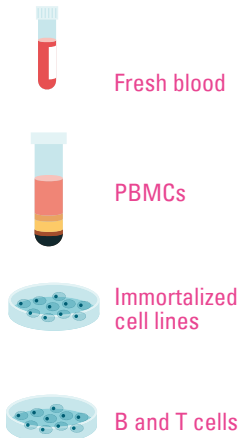

From start to finish

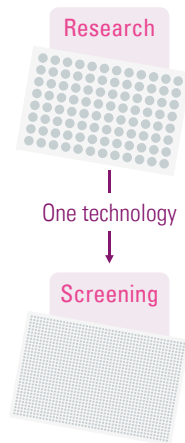

Parallel measurements

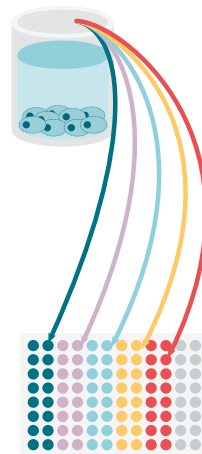


Human G-CSF kit	<i>p.6</i>	Human IL5 kit	<i>p.19</i>	Human & Mouse TGFβ1 kit	<i>p.32</i>
Human GM-CSF kit (CSF2)	<i>p.7</i>	Human IL6 kit	<i>p.20</i>	Human TNFα kit	<i>p.33</i>
Human IFNα kit	<i>p.8</i>	Mouse IL6 kit	<i>p.21</i>	Mouse TNFα kit	<i>p.34</i>
Human IFNβ kit	<i>p.9</i>	Human IL8 kit	<i>p.22</i>	Cynomolgus TNFα kit	<i>p.35</i>
Human IFNγ kit	<i>p.10</i>	Human IL10 kit	<i>p.23</i>	Human CCL2 kit (MCP1)	<i>p.36</i>
Mouse IFNγ kit	<i>p.11</i>	Human IL12 kit	<i>p.24</i>	Mouse CCL2 kit (MCP1)	<i>p.37</i>
Human IL1α kit	<i>p.12</i>	Human IL13 kit	<i>p.25</i>	Human CCL3 kit (MIP1α)	<i>p.38</i>
Mouse IL1α kit	<i>p.13</i>	Human IL17 kit	<i>p.26</i>	Human CCL4 kit (MIP1β)	<i>p.39</i>
Human IL1β kit	<i>p.14</i>	Mouse IL17 kit	<i>p.27</i>	Human CCL5 kit (RANTES)	<i>p.40</i>
Mouse IL1β kit	<i>p.15</i>	Human IL18 kit	<i>p.28</i>	Human CCL20 kit	<i>p.41</i>
Human IL2 kit	<i>p.16</i>	Human IL22 kit	<i>p.29</i>	Human CXCL1 kit (GROα)	<i>p.42</i>
Mouse IL2 kit	<i>p.17</i>	Human IL23 kit	<i>p.30</i>	Human CXCL10 kit (IP10)	<i>p.43</i>
Human IL4 kit	<i>p.18</i>	Human IL31 kit	<i>p.31</i>		

Human G-CSF kit

	Diluent	DMEM	RPMI
Assay range	47 to 4,500 pg/mL		
LOD	27 pg/mL	27 pg/mL	27 pg/mL
LOQ	47 pg/mL	-	-
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (09/136) value (IU/mL) = 0,07 x HTRF hGCSF value (pg/mL)		

Ordering information:

Product	Size	Part#
Human G-CSF kit	500 tests	62HGCSFPEG
Human G-CSF kit	10,000 tests	62HGCSFPEH

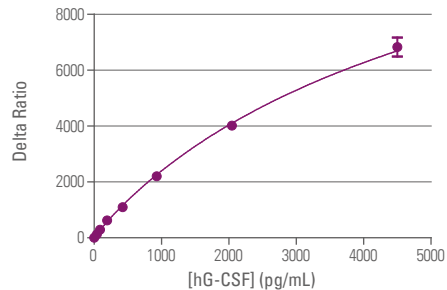


Human GM-CSF kit (CSF2)

	Diluent	DMEM	RPMI
Assay range	54 to 8,500 pg/mL		
LOD	29 pg/mL	29 pg/mL	26 pg/mL
LOQ	54 pg/mL	-	-
Incubation time	3h at room temperature		
Conversion factor	NIBSC (88/646) value (IU/mL) = 0,01 x HTRF hGM-CSF value (pg/mL)		

Ordering information:

Product	Size	Part#
Human GM-CSF kit	500 tests	62HGMCSFPEG
Human GM-CSF kit	10,000 tests	62HGMCSFPEH

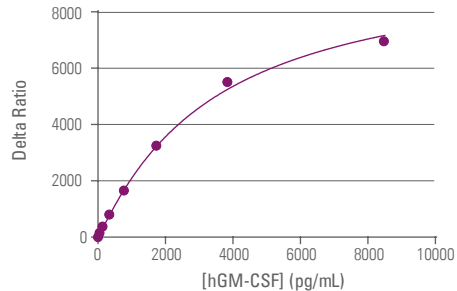


Human IFN α kit

	Diluent	DMEM	RPMI
Assay range	36 to 4000 pg/mL		
LOD	11 pg/mL	24 pg/mL	24 pg/mL
LOQ	36 pg/mL	-	-
Incubation time	2h at room temperature		
Conversion factor	NIBSC (94/784) value (IU/mL) = 0,9 x HTRF hIFN α value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IFN α kit	500 tests	62HIFNAPEG
Human IFN α kit	10,000 tests	62HIFNAPEH

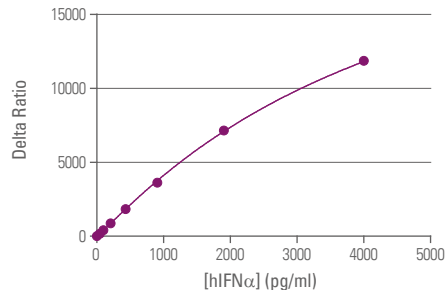


Human IFN β kit

	Diluent	DMEM	RPMI
Assay range	19 to 4,000 pg/mL		
LOD	7 pg/mL	21 pg/mL	20 pg/mL
LOQ	19 pg/mL	-	-
Incubation time	3h at room temperature		
Conversion factor	NIBSC (00/572) value (IU/mL) = 0,44 x HTRF hIFN β value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IFN β kit	500 tests	62HIFNBPEG
Human IFN β kit	10,000 tests	62HIFNBPEH

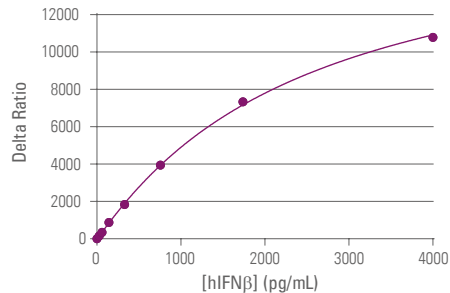


Human IFN γ kit

	Diluent	DMEM	RPMI
Assay range	21 to 4,000 pg/mL		
LOD	14 pg/mL	30 pg/mL	34 pg/mL
LOQ	21 pg/mL	-	-
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (82/587) value (IU/mL) = 0,019 x HTRF hIFN γ value (pg/mL)		


Ordering information:

Product	Size	Part#
Human IFN γ kit	1 x 96 tests	62HIFNGPET
Human IFN γ kit	500 tests	62HIFNGPEG
Human IFN γ kit	10,000 tests	62HIFNGPEH

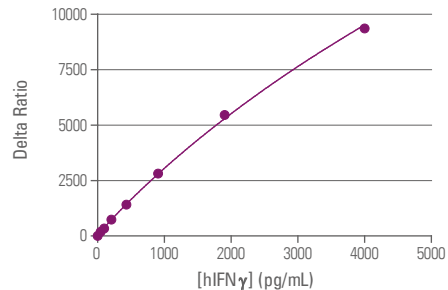

Mouse IFN γ kit

	Diluent	DMEM	RPMI
Assay range	34 to 3,000 pg/mL		
LOD	23 pg/mL	31 pg/mL	30 pg/mL
LOQ	34 pg/mL	-	-
Incubation time	2h at room temperature		

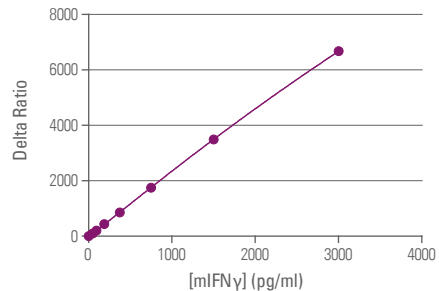

Ordering information:


Product	Size	Part#
Mouse IFN γ kit	500 tests	62MIFNGPEG
Mouse IFN γ kit	10,000 tests	62MIFNGPEH

Human IL1 α kit

	Diluent	DMEM	RPMI
Assay range	20.5 to 5,000 pg/mL		
LOD	8 pg/mL	7 pg/mL	7 pg/mL
LOQ	20.5 pg/mL	-	-
Incubation time	2h at room temperature		
Conversion factor	NIBSC (86/632) value (IU/mL) = 10 x HTRF hIL1 α value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL1 α kit	500 tests	62HIL1APEG
Human IL1 α kit	10,000 tests	62HIL1APEH

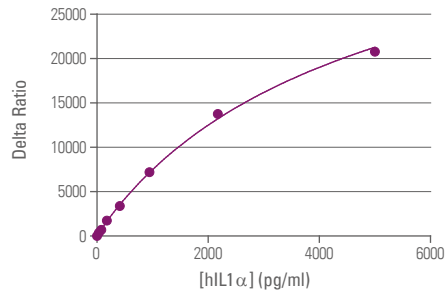


Mouse IL1 α kit

	Diluent	DMEM	RPMI
Assay range	62 to 4,000 pg/mL		
LOD	34 pg/mL	37 pg/mL	41 pg/mL
LOQ	62 pg/mL	-	-
Incubation time	3h at room temperature		
Conversion factor	NIBSC (93/672) value (IU/mL) = 1,2 x HTRF mL1 α value (pg/mL)		

Ordering information:

Product	Size	Part#
Mouse IL1 α kit	500 tests	62MIL1APEG
Mouse IL1 α kit	10,000 tests	62MIL1APEH

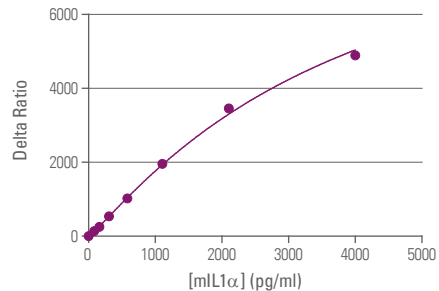


Human IL1 β kit

	Diluent	DMEM	RPMI
Assay range	39 to 6,500 pg/mL		
LOD	5 pg/mL	21 pg/mL	6 pg/mL
LOQ	39 pg/mL	-	-
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (86/680) value (IU/mL) = 0,1 x HTRF hIL1 β value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL1 β kit	1 x 96 tests	62HIL1BPET
Human IL1 β kit	500 tests	62HIL1BPEG
Human IL1 β kit	10,000 tests	62HIL1BPEH

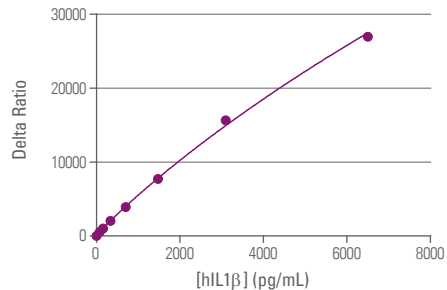


Mouse IL1 β kit

	Diluent	DMEM	RPMI
Assay range	46 to 4,200 pg/mL		
LOD	14 pg/mL	39 pg/mL	37 pg/mL
LOQ	46 pg/mL	-	-
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (93/668) value (IU/mL) = 0,1 x HTRF mL1 β value (pg/mL)		

Ordering information:

Product	Size	Part#
Mouse IL1 β kit	500 tests	62MIL1BPEG
Mouse IL1 β kit	10,000 tests	62MIL1BPEH

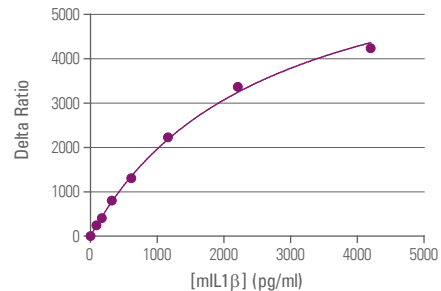


Human IL2 kit

	Diluent	DMEM	RPMI
Assay range	37 to 8,000 pg/mL		
LOD	7 pg/mL	27 pg/mL	17 pg/mL
LOQ	37 pg/mL	-	-
Incubation time	3h at room temperature		
Conversion factor	NIBSC (86/500) value (IU/mL) = 0,01 x HTRF hIL2 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL2 kit	1 x 96 tests	62HIL02PET
Human IL2 kit	500 tests	62HIL02PEG
Human IL2 kit	10,000 tests	62HIL02PEH

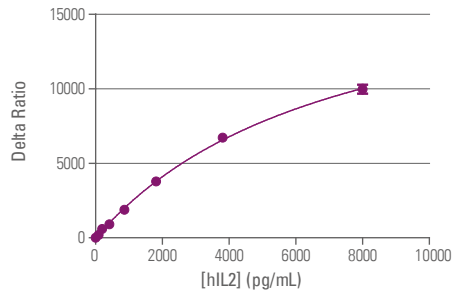


Mouse IL2 kit

	Diluent	DMEM	RPMI
Assay range	35 to 2,560 pg/mL		
LOD	12 pg/mL	27 pg/mL	16 pg/mL
LOQ	35 pg/mL	-	-
Incubation time	Overnight at room temperature		

Ordering information:

Product	Size	Part#
Mouse IL2 kit	500 tests	62MIL02PEG
Mouse IL2 kit	10,000 tests	62MIL02PEH

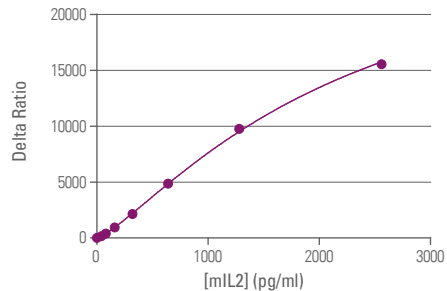


Human IL4 kit

	Diluent	DMEM	RPMI
Assay range	29 to 4,000 pg/mL		
LOD	8 pg/mL	19 pg/mL	10 pg/mL
LOQ	29 pg/mL	-	-
Incubation time	2h at room temperature		
Conversion factor	NIBSC (88/656) value (IU/mL) = 0,012 x HTRF hIL4 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL4 kit	500 tests	62HIL04PEG
Human IL4 kit	10,000 tests	62HIL04PEH

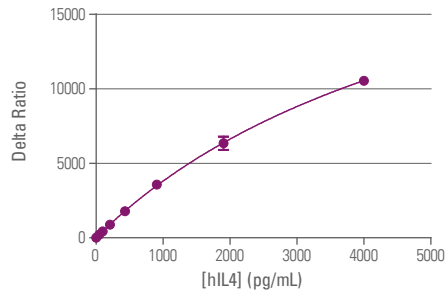


Human IL5 kit

	Diluent	DMEM	RPMI
Assay range	71 to 10,000 pg/mL		
LOD	15 pg/mL	24 pg/mL	14 pg/mL
LOQ	71 pg/mL	-	-
Incubation time	2h at room temperature		
Conversion factor	NIBSC (90/586) value (IU/mL) = 0,012 x HTRF hIL5 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL5 kit	500 tests	62HIL05PEG
Human IL5 kit	10,000 tests	62HIL05PEH

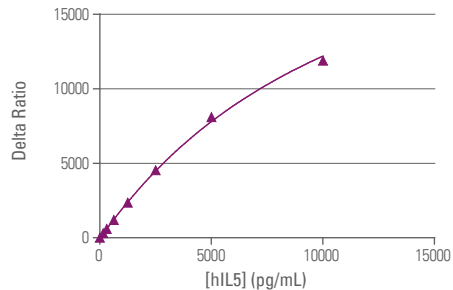


Human IL6 kit

	Diluent	DMEM	RPMI
Assay range	34 to 7,500 pg/mL		
LOD	8 pg/mL	19 pg/mL	23 pg/mL
LOQ	34 pg/mL	-	-
Incubation time	2h at room temperature		
Conversion factor	NIBSC (89/548) value (IU/mL) = 0,1 x HTRF hIL6 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL6 kit	1 x 96 tests	62HIL06PET
Human IL6 kit	500 tests	62HIL06PEG
Human IL6 kit	10,000 tests	62HIL06PEH

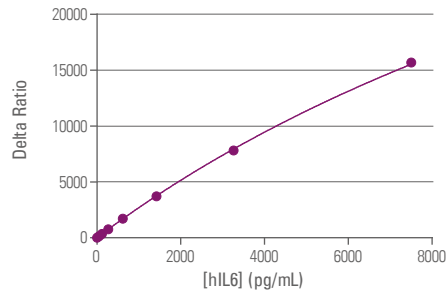


Mouse IL6 kit

	Diluent	DMEM	RPMI
Assay range	19 to 7,000 pg/mL		
LOD	6 pg/mL	13 pg/mL	9 pg/mL
LOQ	19 pg/mL	-	-
Incubation time	1h at room temperature		
Conversion factor	NIBSC (93/730) value (IU/mL) = 0,7 x HTRF mL6 value (pg/mL)		

Ordering information:

Product	Size	Part#
Mouse IL6 kit	500 tests	62MIL06PEG
Mouse IL6 kit	10,000 tests	62MIL06PEH

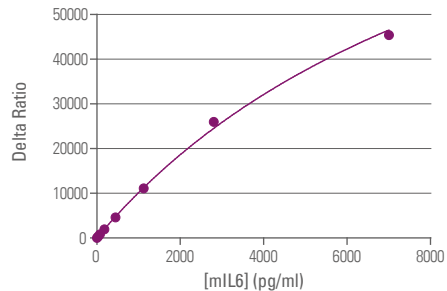


Human IL8 kit

	Diluent	DMEM	RPMI
Assay range	32 to 4,000 pg/mL		
LOD	6 pg/mL	11 pg/mL	8 pg/mL
LOQ	32 pg/mL	-	-
Incubation time	1h at room temperature		
Conversion factor	NIBSC (89/520) value (IU/mL) = 0,01 x HTRF hIL8 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL8 kit	1 x 96 tests	62HIL08PET
Human IL8 kit	500 tests	62HIL08PEG
Human IL8 kit	10,000 tests	62HIL08PEH

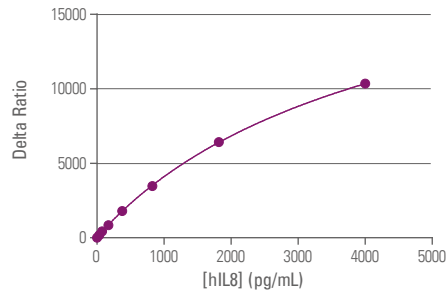


Human IL10 kit

	Diluent	DMEM	RPMI
Assay range	22 to 6,000 pg/mL		
LOD	7 pg/mL	13 pg/mL	8 pg/mL
LOQ	22 pg/mL	-	-
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (93/722) value (IU/mL) = 0,005 x HTRF hIL10 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL10 kit	500 tests	62HIL10PEG
Human IL10 kit	10,000 tests	62HIL10PEH

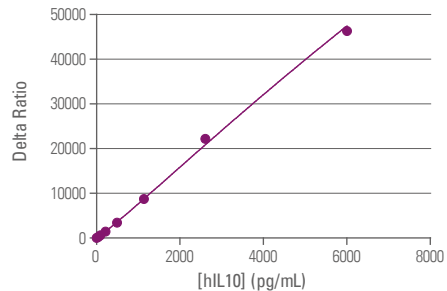


Human IL12 kit

	Diluent	DMEM	RPMI
Assay range	81 to 13,000 pg/mL		
LOD	14 pg/mL	61 pg/mL	34 pg/mL
LOQ	81 pg/mL	-	-
Incubation time	2h at room temperature		
Conversion factor	NIBSC (95/544) value (IU/mL) = 0,01 x HTRF hIL12 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL12 kit	500 tests	62HIL12PEG
Human IL12 kit	10,000 tests	62HIL12PEH

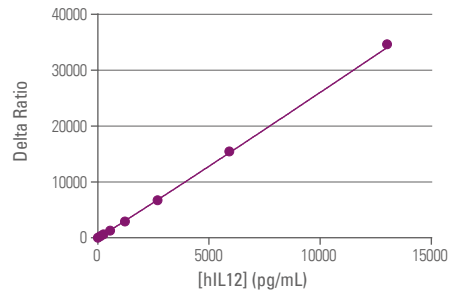


Human IL13 kit

	Diluent	DMEM	RPMI
Assay range	17 to 5,000 pg/mL		
LOD	6 pg/mL	19 pg/mL	23 pg/mL
LOQ	17 pg/mL	24 pg/mL	25 pg/mL
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (94/622) value (IU/mL) = $1.2 \cdot 10^{-3} \times$ HTRF hIL13 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL13 kit	500 tests	62HIL13PEG
Human IL13 kit	10,000 tests	62HIL13PEH

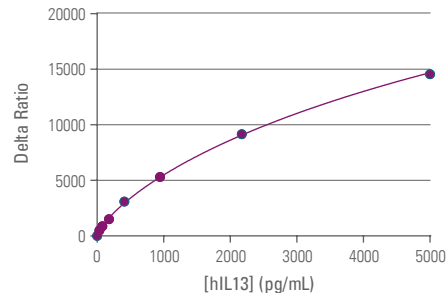


Human IL17 kit

	Diluent	DMEM	RPMI
Assay range	33 to 10,000 pg/mL		
LOD	7 pg/mL	22 pg/mL	22 pg/mL
LOQ	33 pg/mL	-	-
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (01/420) value (IU/mL) = 0,02 x HTRF hIL17 value (pg/mL)		


Ordering information:

Product	Size	Part#
Human IL17 kit	1 x 96 tests	62HIL17PET
Human IL17 kit	500 tests	62HIL17PEG
Human IL17 kit	10,000 tests	62HIL17PEH

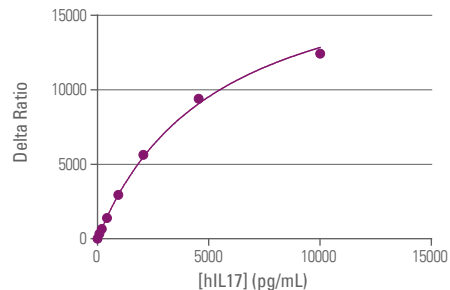

Mouse IL17 kit

	Diluent	DMEM	RPMI
Assay range	16 to 4,000 pg/mL		
LOD	4 pg/mL	9 pg/mL	6 pg/mL
LOQ	16 pg/mL	-	-
Incubation time	1h at room temperature		

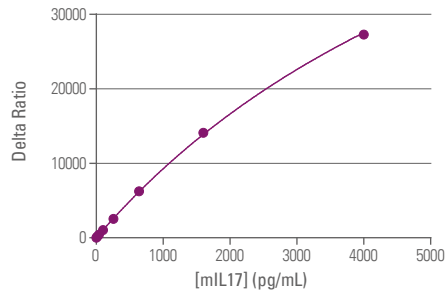

Ordering information:


Product	Size	Part#
Mouse IL17 kit	500 tests	62MIL17PEG
Mouse IL17 kit	10,000 tests	62MIL17PEH

Human IL18 kit

	Diluent	DMEM	RPMI
Assay range	43 to 8,000 pg/mL		
LOD	18 pg/mL	32 pg/mL	84 pg/mL
LOQ	43 pg/mL	-	-
Incubation time	1h at room temperature		
Conversion factor	NIBSC (03/200) value (IU/mL) = 5 x HTRF hIL18 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human IL18 kit	500 tests	62HIL18PEG
Human IL18 kit	10,000 tests	62HIL18PEH

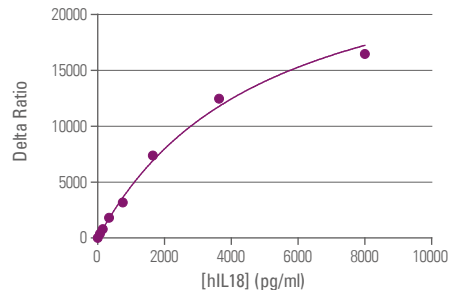


Human IL22 kit

	Diluent	DMEM	RPMI
Assay range	26 to 5,000 pg/mL		
LOD	9 pg/mL	16 pg/mL	13 pg/mL
LOQ	26 pg/mL	35 pg/mL	28 pg/mL
Incubation time	Overnight at room temperature (reading possible from 3h)		
Conversion factor	No International standard available		


Ordering information:

Product	Size	Part#
Human IL22 kit	500 tests	62HIL22PEG
Human IL22 kit	10,000 tests	62HIL22PEH

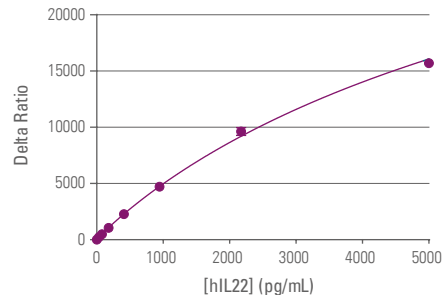

Human IL23 kit

	Diluent	DMEM	RPMI
Assay range	92 to 30,000 pg/mL		
LOD	30 pg/mL	46 pg/mL	40 pg/mL
LOQ	92 pg/mL	-	-
Incubation time	3h at 4°C		

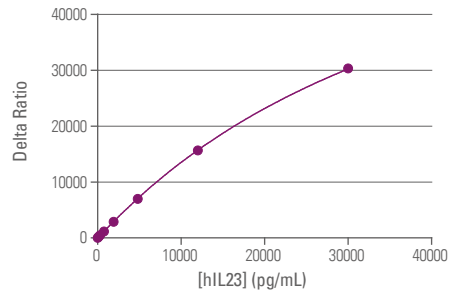


Ordering information:

Product	Size	Part#
Human IL23 kit	500 tests	62HIL23PEG
Human IL23 kit	10,000 tests	62HIL23PEH

Human IL31 kit

	Diluent	DMEM	RPMI
Assay range	87 to 20,000 pg/mL		
LOD	18 pg/mL	50 pg/mL	36 pg/mL
LOQ	87 pg/mL	-	-
Incubation time	Overnight at room temperature		

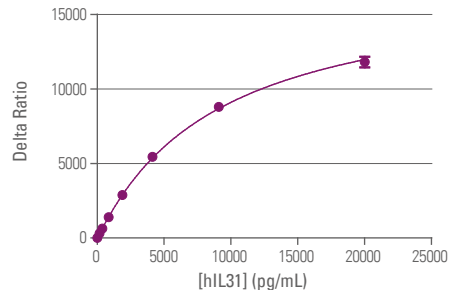

Ordering information:


Product	Size	Part#
Human IL31 kit	500 tests	62HIL31PEG
Human IL31 kit	10,000 tests	62HIL31PEH

Human & Mouse TGFβ1 kit

	Diluent	DMEM + 4% FCS	RPMI + 4% FCS
Assay range	19 to 2,000 pg/mL		
LOD	4 pg/mL	13 pg/mL	6 pg/mL
LOQ	19 pg/mL	26 pg/mL	28 pg/mL
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (89/514) value (U/mL) = 0,017 x HTRF hTGFβ1 value (pg/mL)		

Ordering information:

Product	Size	Part#
TGFβ1 kit	500 tests	62HTGFBPEG
TGFβ1 kit	10,000 tests	62HTGFBPEH

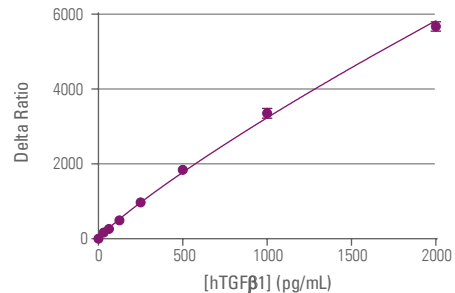


Human TNF α kit

	Diluent	DMEM	RPMI
Assay range	29 to 2,500 pg/mL		
LOD	5 pg/mL	17 pg/mL	11 pg/mL
LOQ	29 pg/mL	-	-
Incubation time	2h at room temperature		
Conversion factor	NIBSC (12/154) value (IU/mL) = 0,1 x HTRF hTNF α value (pg/mL)		


Ordering information:

Product	Size	Part#
Human TNF α kit	1 x 96 tests	62HTNFAPET
Human TNF α kit	500 tests	62HTNFAPEG
Human TNF α kit	10,000 tests	62HTNFAPEH

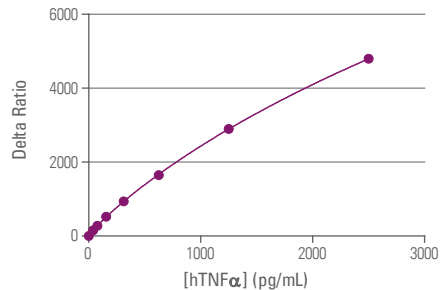

Mouse TNF α kit

	Diluent	DMEM	RPMI
Assay range	20 to 6,000 pg/mL		
LOD	3 pg/mL	11 pg/mL	8 pg/mL
LOQ	20 pg/mL	-	-
Incubation time	1h at room temperature		

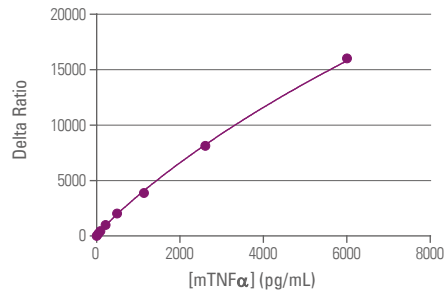


Ordering information:

Product	Size	Part#
Mouse TNF α kit	500 tests	62MTNFAPEG
Mouse TNF α kit	10,000 tests	62MTNFAPEH

Cynomolgus TNF α kit

	Diluent	DMEM	RPMI
Assay range	27 to 3,000 pg/mL		
LOD	12 pg/mL	44 pg/mL	40 pg/mL
LOQ	27 pg/mL	-	-
Incubation time	Overnight at room temperature		

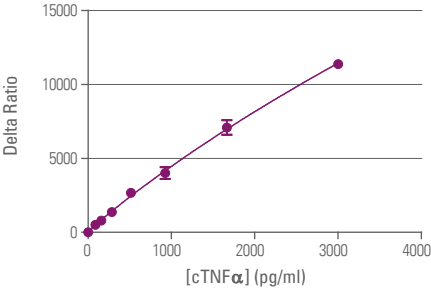

Ordering information:


Product	Size	Part#
Cynomolgus TNF α kit	500 tests	62CTNFAPEG
Cynomolgus TNF α kit	10,000 tests	62CTNFAPEH

Human CCL2 kit (MCP1)

	Diluent	DMEM	RPMI
Assay range	90 to 3,000 pg/mL		
LOD	20 pg/mL	28 pg/mL	20 pg/mL
LOQ	90 pg/mL	-	-
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (92/794) value (IU/mL) = 0,9 x HTRF hCCL2 value (pg/mL)		


Ordering information:

Product	Size	Part#
Human CCL2 kit (MCP1)	500 tests	62HCCL2PEG
Human CCL2 kit (MCP1)	10,000 tests	62HCCL2PEH

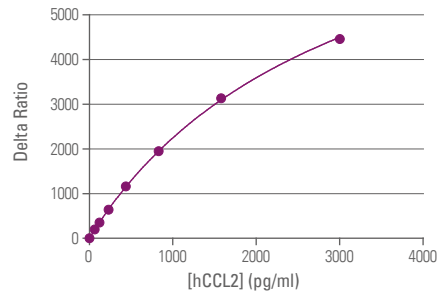

Mouse CCL2 kit (MCP1)

	Diluent	DMEM	RPMI
Assay range	13 to 1,800 pg/mL		
LOD	6 pg/mL	18 pg/mL	17 pg/mL
LOQ	13 pg/mL	-	-
Incubation time	Overnight at room temperature		

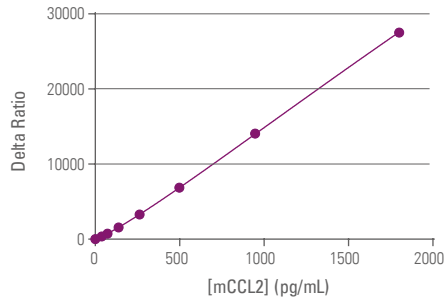

Ordering information:


Product	Size	Part#
Mouse CCL2 kit	500 tests	62MCCL2PEG
Mouse CCL2 kit	10,000 tests	62MCCL2PEH

Human CCL3 kit (MIP1 α)

	Diluent	DMEM	RPMI
Assay range	58 to 4,400 pg/mL		
LOD	28 pg/mL	35 pg/mL	25 pg/mL
LOQ	58 pg/mL	52 pg/mL	56 pg/mL
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (92/518) value (IU/mL) = 0,00005 x HTRF hCCL3 value (pg/mL)		

Ordering information:

Product	Size	Part#
Human CCL3 kit	500 tests	62HCCL3PEG
Human CCL3 kit	10,000 tests	62HCCL3PEH

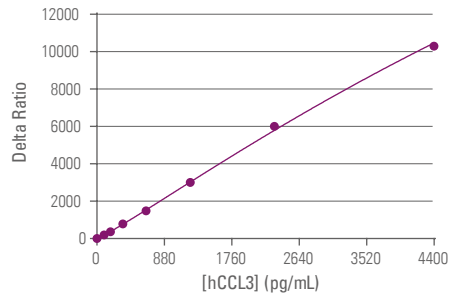


Human CCL4 kit (MIP1 β)

	Diluent	DMEM	RPMI
Assay range	20 to 5,000 pg/mL		
LOD	7 pg/mL	9 pg/mL	10 pg/mL
LOQ	20 pg/mL	-	-
Incubation time	1h at room temperature		

Ordering information:

Product	Size	Part#
Human CCL4 kit (MIP1 β)	500 tests	62HCCL4PEG
Human CCL4 kit (MIP1 β)	10,000 tests	62HCCL4PEH

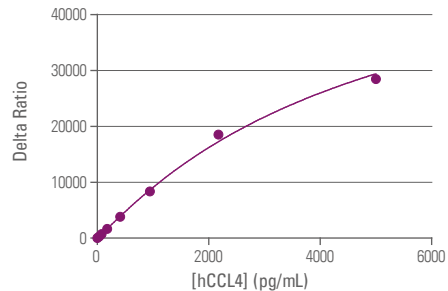


Human CCL5 kit (RANTES)

	Diluent	DMEM	RPMI
Assay range	18 to 2,000 pg/mL		
LOD	7 pg/mL	12 pg/mL	9 pg/mL
LOQ	18 pg/mL	-	-
Incubation time	Overnight at room temperature		
Conversion factor	NIBSC (92/520) value (IU/mL) = 0,015 x HTRF hCCL5 value (pg/mL)		


Ordering information:

Product	Size	Part#
Human CCL5 kit	500 tests	62HCCL5PEG
Human CCL5 kit	10,000 tests	62HCCL5PEH

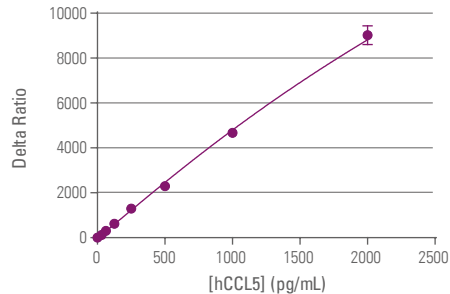

Human CCL20 kit

	Diluent	DMEM	RPMI
Assay range	47 to 2,000 pg/mL		
LOD	15 pg/mL	82 pg/mL	34 pg/mL
LOQ	47 pg/mL	-	-
Incubation time	Overnight at room temperature		

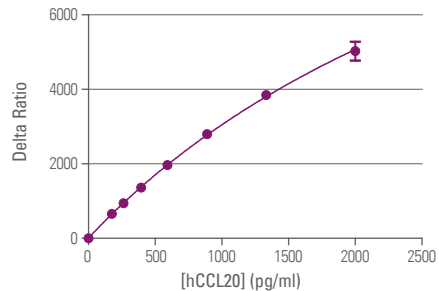


Ordering information:

Product	Size	Part#
Human CCL20 kit	500 tests	62HCC20PEG
Human CCL20 kit	10,000 tests	62HCC20PEH

Human CXCL1 kit (GRO α)

	Diluent	DMEM	RPMI
Assay range	36 to 4,000 pg/mL		
LOD	13 pg/mL	33 pg/mL	40 pg/mL
LOQ	36 pg/mL	-	-
Incubation time	2h at 4°C		

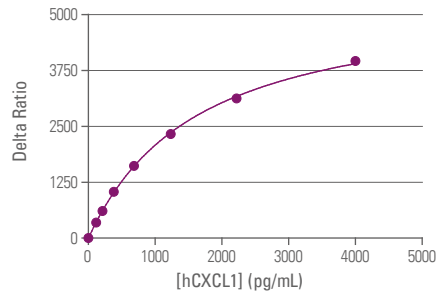

Ordering information:


Product	Size	Part#
Human CXCL1 kit (GRO α)	500 tests	62HCXC1PEG
Human CXCL1 kit (GRO α)	10,000 tests	62HCXC1PEH

Human CXCL10 kit (IP10)

	Diluent	DMEM	RPMI
Assay range	26 to 5,000 pg/mL		
LOD	1.8 pg/mL	3.8 pg/mL	4.2 pg/mL
LOQ	7 pg/mL	-	-
Incubation time	2h at room temperature		

Ordering information:

Product	Size	Part#
Human CXCL10 kit (IP10)	500 tests	62HCX10PEG
Human CXCL10 kit (IP10)	10,000 tests	62HCX10PEH

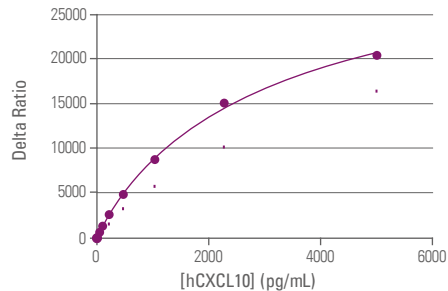

MORE THAN PRODUCTS, KNOWLEDGE!

Find helpful reference material, content, and expert advice by visiting our website at www.cisbio.com/cytokines





VIDEOS

| How to quantify secreted cytokines using no-wash, add-and-read HTRF assays

| How to combine Cisbio assays to measure several cytokines from the same supernatant



APPLICATION NOTES

| Guidelines from PBMC isolation to cytokine assay optimisation

| HTRF® measurement of cytokine release from fresh blood samples

POSTER

| Cytokine secretion profile on PBMCs



Available assays listed by species

Human G-CSF kit	<i>p.6</i>
Human GM-CSF kit (CSF2)	<i>p.7</i>
Human IFN α kit	<i>p.8</i>
Human IFN β kit	<i>p.9</i>
Human IFN γ kit	<i>p.10</i>
Human IL1 α kit	<i>p.12</i>
Human IL1 β kit	<i>p.14</i>
Human IL2 kit	<i>p.16</i>
Human IL4 kit	<i>p.18</i>
Human IL5 kit	<i>p.19</i>
Human IL6 kit	<i>p.20</i>
Human IL8 kit	<i>p.22</i>
Human IL10 kit	<i>p.23</i>
Human IL12 kit	<i>p.24</i>
Human IL13 kit	<i>p.25</i>

Human IL17 kit	<i>p.26</i>
Human IL18 kit	<i>p.28</i>
Human IL22 kit	<i>p.29</i>
Human IL23 kit	<i>p.30</i>
Human IL31 kit	<i>p.31</i>
Human & Mouse TGF β 1 kit	<i>p.32</i>
Human TNF α kit	<i>p.33</i>
Human CCL2 kit (MCP1)	<i>p.36</i>
Human CCL3 kit (MIP1 α)	<i>p.38</i>
Human CCL4 kit (MIP1 β)	<i>p.39</i>
Human CCL5 kit (RANTES)	<i>p.40</i>
Human CCL20 kit	<i>p.41</i>
Human CXCL1 kit (GRO α)	<i>p.42</i>
Human CXCL10 kit (IP10)	<i>p.43</i>

Mouse IFN γ kit	<i>p.11</i>
Mouse IL1 α kit	<i>p.13</i>
Mouse IL1 β kit	<i>p.15</i>
Mouse IL2 kit	<i>p.17</i>
Mouse IL6 kit	<i>p.21</i>
Mouse IL17 kit	<i>p.27</i>
Human & Mouse TGF β 1 kit	<i>p.32</i>
Mouse TNF α kit	<i>p.34</i>
Mouse CCL2 kit (MCP1)	<i>p.37</i>

Cynomolgus TNF α kit	<i>p.35</i>
-----------------------------	-------------

Europe and other countries
+33(0)466 796 705 | info@cisbio.com

U.S. and Canada
1-888-963-4567 | htrfinfo@cisbio.us

China
+86 21 5018 9880 | htrfinfoasia@cisbio.cn

Japan
+81-(0)43-306-8712 | htrf@cisbio.jp

www.cisbio.com



cisbio